

Automated set-up of the Lonza PyroGene[®] Recombinant Factor C Assay Endotoxin Detection Test on the ASSIST PLUS

Introduction

The Lonza PyroGene[®] Recombinant Factor C (rFC) Endotoxin Detection Assay is an alternative to the traditional limulus amoebocyte lysate (LAL) assay, which is widely used to screen for bacterial endotoxin contamination in human and animal parenteral pharmaceuticals and medical devices. The rFC test is used in both high and low throughput laboratories and, unlike the LAL assay, it is not derived from horseshoe crab blood. Setting up the test requires manual preparation of 10-fold diluted standards from the endotoxin stock solution supplied. Standards and samples are added in duplicate wells to a 96 well plate. To check for product inhibition, positive product controls (PPCs), which are samples spiked with a known concentration of endotoxin, are tested alongside samples. Following the initial plating of the standards, samples and PPCs, a 10-minute pre-incubation is performed,


during which the user prepares a working solution that consists of the fluorogenic substrate, assay buffer and rFC enzyme.

This application note demonstrates that the preparation of the standards, samples, PPC samples and blanks, and their addition to the 96 well plate, can be easily automated on the ASSIST PLUS pipetting robot using a D-ONE single channel pipetting module. Addition of the working reagent is then completed using the VOYAGER adjustable tip spacing multichannel pipette. Automation of all the pipetting steps reduces the opportunity for pipetting errors, and ensures robust reproducibility. The key quality indicators in this assay are the correlation coefficient of the standard curve and coefficient of sample variation (CV).

Key benefits:

- Automated preparation of standard dilutions eliminates pipetting errors that would invalidate entire runs.
- Users can perform testing of full plates or partial plates by using the D-ONE single channel pipetting module in combination with the flexible and easy to program ASSIST PLUS.
- Error-free pipetting ensures replicate samples with tight CV values, reducing the likelihood of repeat testing.
- On-screen prompts guide the user through instrument set-up.

Overview: How to perform the PyroGene[®] rFC endotoxin detection assay

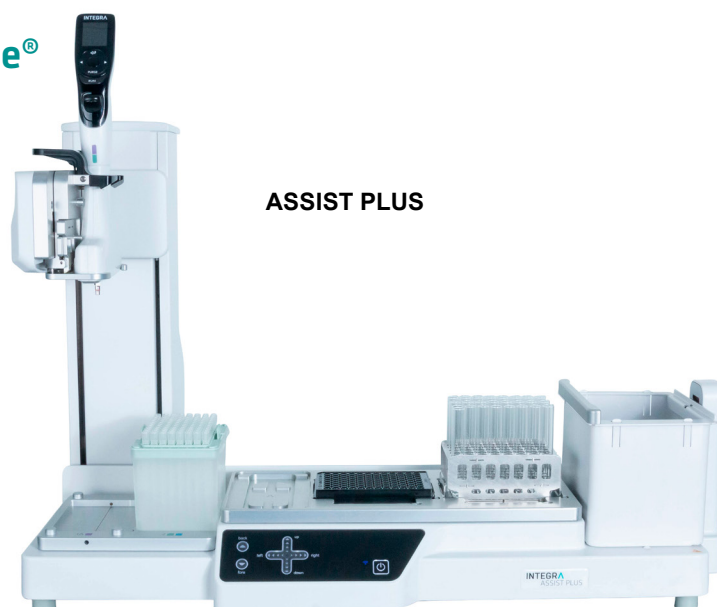

The PyroGene® Recombinant Factor C (rFC) Endotoxin Detection Assay is a fluorogenic assay that requires a fluorescence microplate reader – such as PyroWave® XM Reader paired with WinKQCL® Endotoxin Detection & Analysis Software – to measure endotoxin values. Prior to setting up the plate, a template is prepared in WinKQCL® Software. **Figure 1** shows the WinKQCL® Software template that was developed for use with samples plated by the ASSIST PLUS. This template allows up to 21 samples to be tested in duplicate on one plate with paired PPC samples. It is designed to provide the most flexibility for varying numbers of samples, while still allowing use of an 8 channel multichannel pipette to deliver rFC working solution to the plate, as written in the instructions for use.

	1	2	3	4	5	6	7	8	9	10	11	12
A	BLNK	BLNK	543 Filler 230301	543 Filler 230301	S11 S6 Test .25 230301	S11 S6 Test .25 230301	S12+ S6 Test .25 230301	S12+ S6 Test .25 230301	S27 S14 Test .25 230301	S27 S14 Test .25 230301	S28+ S14 Test .25 230301	S28+ S14 Test .25 230301
B	STD1 0.005	STD1 0.005	STD2 0.05	STD2 0.05	S13 S7 Test .025 230301	S13 S7 Test .025 230301	S14+ S7 Test .025 230301	S14+ S7 Test .025 230301	S29 S15 Test .025 230301	S29 S15 Test .025 230301	S30+ S15 Test .025 230301	S30+ S15 Test .025 230301
C	STD3 0.5	STD3 0.5	STD4 5	STD4 5	S15 S8 Test 0 230301	S15 S8 Test 0 230301	S16+ S8 Test 0 230301	S16+ S8 Test 0 230301	S31 S16 Test 0 230301	S31 S16 Test 0 230301	S32+ S16 Test 0 230301	S32+ S16 Test 0 230301
D	S1 S1 Test 2.5 230301	S1 S1 Test 2.5 230301	S2+ S1 Test 2.5 230301	S2+ S1 Test 2.5 230301	S17 S9 Test 2.5 230301	S17 S9 Test 2.5 230301	S18+ S9 Test 2.5 230301	S18+ S9 Test 2.5 230301	S33 S17 Test 2.5 230301	S33 S17 Test 2.5 230301	S34+ S17 Test 2.5 230301	S34+ S17 Test 2.5 230301
E	S3 S2 Test .25 230301	S3 S2 Test .25 230301	S4+ S2 Test .25 230301	S4+ S2 Test .25 230301	S19 S10 Test .25 230301	S19 S10 Test .25 230301	S20+ S10 Test .25 230301	S20+ S10 Test .25 230301	S35 S18 Test .25 230301	S35 S18 Test .25 230301	S36+ S18 Test .25 230301	S36+ S18 Test .25 230301
F	S5 S3 Test .025 230301	S5 S3 Test .025 230301	S6+ S3 Test .025 230301	S6+ S3 Test .025 230301	S21 S11 Test .025 230301	S21 S11 Test .025 230301	S22+ S11 Test .025 230301	S22+ S11 Test .025 230301	S37 S19 Test .025 230301	S37 S19 Test .025 230301	S38+ S19 Test .025 230301	S38+ S19 Test .025 230301
G	S7 S4 Test 0 230301	S7 S4 Test 0 230301	S8+ S4 Test 0 230301	S8+ S4 Test 0 230301	S23 S12 Test 0 230301	S23 S12 Test 0 230301	S24+ S12 Test 0 230301	S24+ S12 Test 0 230301	S39 S20 Test 0 230301	S39 S20 Test 0 230301	S40+ S20 Test 0 230301	S40+ S20 Test 0 230301
H	S9 S5 Test 2.5 230301	S9 S5 Test 2.5 230301	S10+ S5 Test 2.5 230301	S10+ S5 Test 2.5 230301	S25 S13 Test 2.5 230301	S25 S13 Test 2.5 230301	S26+ S13 Test 2.5 230301	S26+ S13 Test 2.5 230301	S41 S21 Test 0 230301	S41 S21 Test 0 230301	S42+ S21 Test 0 230301	S42+ S21 Test 0 230301

Figure 1: Lonza rFC template in WinKQCL® Software. Red wells: the position of the standards; Blue wells: samples tested in duplicate; yellow wells: PPC-spiked samples tested in duplicate. Wells A3 and A4 are empty.

The D-ONE 5-300 µl single channel pipetting module is first used to prepare standard dilutions. Following preparation of the standards, PPC is added to the designated wells of the 96 well plate in a repeat dispense mode. Next, standards and samples are added to the plate according to the above template. Standards and samples are added to the plate in duplicate, with samples added to both clean and PPC-spiked wells. Once all samples are added, the plate is pre-incubated at 37 °C for 10 minutes. During the incubation, the user prepares the working reagent in an INTEGRA 10 ml SureFlo™ reagent reservoir. At the end of the 10 minute pre-incubation, the working reagent is added to the plate using a VOYAGER 8 channel 300 µl pipette on the ASSIST PLUS.

TIPS:

- Use pyrogen-free certified GRIPTIPS® in combination with SureFlo reservoirs to ensure accurate results. 10 ml SureFlo reservoirs require a dead volume of less than 30 µl.
- LONG 300 µl GRIPTIPS can access sample volumes of below 1 ml in 13 x 100 mm Lonza pyrogen free test tubes, and will never leak or fall off.
- PPC and samples are dispensed using repeat dispense mode to save time and money.
- Plate layout of the blank, standards and samples in the WinKQCL® Software template is designed to offer the most flexibility for running full or partial plates.

Experimental set-up

Deck position A: LAL water – 10 ml multichannel reagent reservoir

Deck position B: PyroGene® rFC assay plate – 96 well flat bottom plate (Corning)

Deck position C: Standards, samples and PPC spike – INTEGRA tube rack

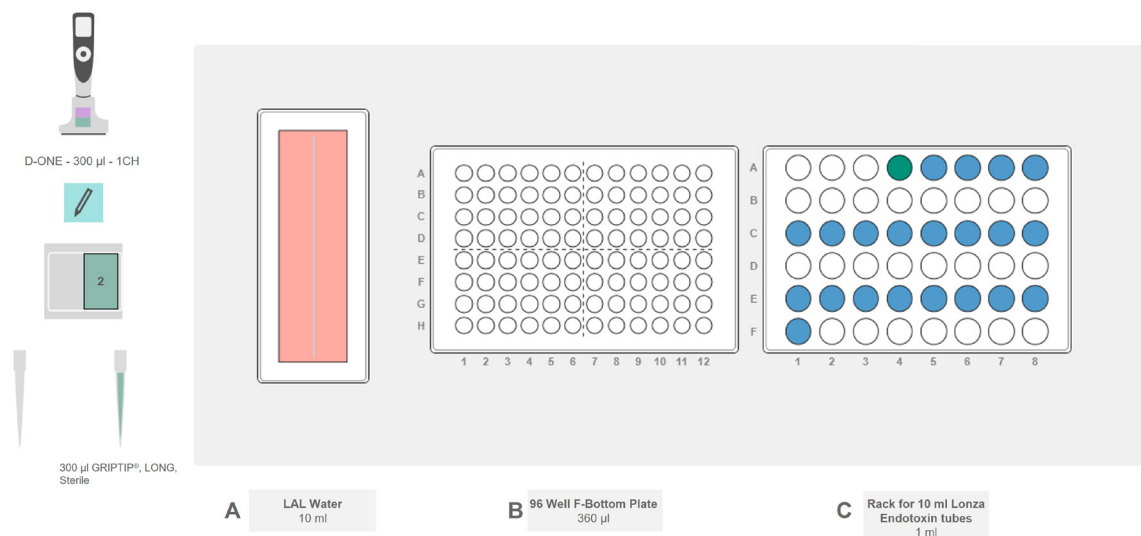


Figure 2: Set-up of the ASSIST PLUS for PyroGene® Assay plate format. **Position A:** LAL water in 10 ml SureFlo reservoir. **Position B:** 96 well assay plate. **Position C:** Tube rack with Lonza pyrogen-free tubes (empty, magenta), 250 µl of 20 EU/ml endotoxin standard (green) and 1 ml of sample (blue).

1. PyroGene® rFC assay plate set-up

STEP: Standards, samples and PPC spiked samples are added to the plate.

HOW TO: Pair the D-ONE 0.5-300 µl single channel pipetting module with the ASSIST PLUS pipetting robot. Place pyrogen-free test tubes containing 1 ml sample in the tube rack on deck position C (blue tubes in **Figure 2**). Place three empty tubes in positions A1-A3 within the rack. These are used to create the standard dilutions (magenta tubes in **Figure 2**). Place a tube holding 250 µl of 20 EU/ml endotoxin standard in A4 in the rack (green tube in **Figure 2**). Place a 10 ml SureFlo reservoir holding LAL water on deck position A, and a 96 well fluorescence plate on deck position B.

When the 'Lonza PyroGene® Assay plate set-up' program is started, the pipette first dispenses 900 µl of LAL water into each of the three empty dilution tubes (**Figure 3**). Next, 750 µl of LAL water is dispensed into the tube holding 250 µl of 20 EU/ml endotoxin standard, creating a 5 EU/ml standard. Following package insert instructions, a message on the pipette instructs the user to vortex the 5 EU/ml standard for 1 minute. The ASSIST PLUS pauses for the user to perform this step, then restarts when the user acknowledges the message. The next standard is created by transfer of 100 µl of the 5 EU/ml standard to the adjacent tube holding 900 µl of LAL water, followed by vortexing. The remaining two standards are similarly created.



Figure 3: LAL water is dispensed by the D-ONE single channel pipetting module.

Once all the standards are created, 10 μ l of the 5 EU/ml standard is added to each well of the plate designated as a PPC (**Figure 4**). This serves as a control to monitor for sample inhibition of endotoxin detection. Each blank, standard and sample is then added to the appropriate wells in duplicate. Duplicate samples are also added to the PPC wells, creating the 0.5 EU/ml PPC-spiked samples. When all standards and samples have been added to the plate, the plate is pre-incubated at 37 °C for 10 minutes.



Figure 4: PPC is added by the D-ONE single channel pipetting module.

2. PyroGene® assay reagent addition

STEP: Add 100 µl working reagent to each well of the plate.

HOW TO: While the plate is pre-incubating, manually prepare the working reagent in a 10 ml SureFlo reservoir by combining fluorogenic substrate, rFC assay buffer and rFC enzyme solution in a 5:4:1 ratio. Place the working reagent in a clean 10 ml SureFlo reservoir on deck position A (**Figure 5**). Pair the 300 µl 8 channel VOYAGER pipette with the ASSIST PLUS, and exchange the D-ONE tip deck for the standard tip deck. At the end of the 10 minute pre-incubation, place the plate on deck position B. Initiate the VIALAB program 'Lonza PyroGene® Assay reagent addition' to add 100 µl reagent to each well of the plate. When reagent addition is complete, place the plate in the fluorescence microplate reader to complete the assay.

TIPS:

- VIALAB programs can be adapted to accommodate different numbers of samples, giving laboratories total flexibility to meet current and future testing demands.
- LONG 300 µl GRIPTIPS are used in all steps of this assay, simplifying the purchasing process for customers.

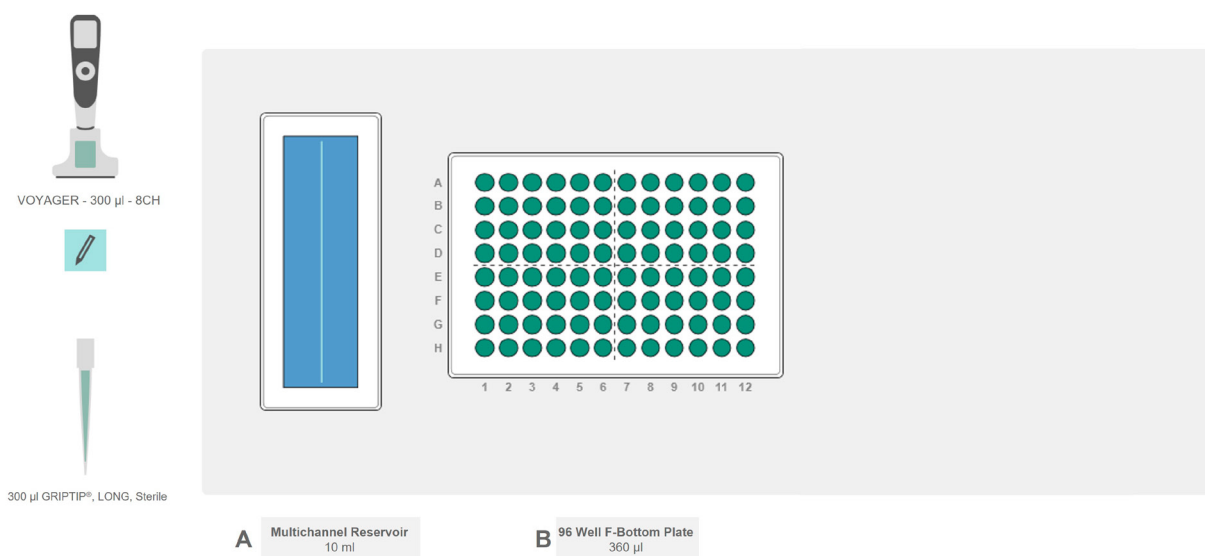


Figure 5: Deck set-up of the ASSIST PLUS for reagent addition. **Position A:** working endotoxin reagent in a 10 ml SureFlo reservoir. **Position B:** 96 well plate containing pre-incubated standards and samples.

Assay verification

Three runs of 21 samples and standards were set up on the ASSIST PLUS and tested according to the PyroGene® rFC Assay instructions for use. Samples consisted of LAL water spiked with known concentrations of endotoxin standard. 5 samples at each concentration were run on each plate, except for the 0 EU/ml sample, which was run 6 times per plate. Data analysis was performed using WinkQCL® Software. **Table 1** shows the concentrations of standards and samples.

Table 1: Endotoxin concentrations of standards and spiked samples.

Standards [EU/ml]	Samples of LAL water spiked with endotoxin [EU/ml]
5	2.5
0.5	0.25
0.05	0.025
0.005	0

Results

Results are displayed in **Tables 2-4**. The standard curve for all runs displayed linearity, and all curves were within the quality parameters as defined by the instructions for use (**Table 2**). Samples spiked with each concentration of endotoxin were detected (**Figure 6**). All unspiked samples remained undetectable at <0.005 EU/ml, which is the cut-off for acceptable endotoxin concentrations in pharmaceuticals and medical devices. All replicates of standards, samples and PPC samples displayed a CV within the acceptable limit of below 25 % (**Table 3**).

Table 2: Standard curve results with quality specifications for runs 1-3.

	Correlation coefficient (0.980-1.000)	Slope (0.760-1.110)	Y-intercept (2.500-5.000)
Run 1	1.000	0.930	4.112
Run 2	0.999	0.903	4.258
Run 3	0.998	0.909	4.267

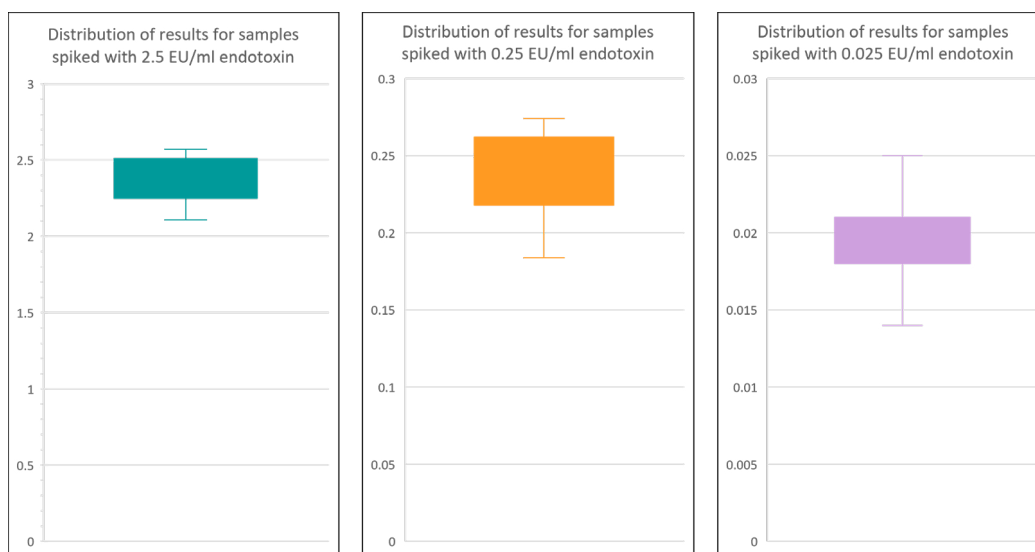


Figure 6: Distribution of results for samples spiked at endotoxin concentrations of 2.5, 0.25 and 0.025 EU/ml.

Table 3: Mean coefficient of variation per plate for paired samples and PPC samples.

	Mean paired sample %CV	Mean PPC %CV	QC fail %CV per plate
Run 1	4.19	4.37	0
Run 2	1.84	4.03	0
Run 3	3.66	3.87	0
Overall Mean	3.23	4.09	0

Remarks

- **Partial plates:** The supplied VIALAB programs can be adapted for partial plates, or for running samples in triplicate.
- **Run report:** If the ASSIST PLUS pipetting robot is connected to the PC, VIALAB programs can be started directly from the PC. After the run, a report is automatically generated, documenting details such as start and end times, user identification, calculated volumes and any errors that occurred. This offers a convenient way to fulfill regulatory requirements.

Conclusion

- The ASSIST PLUS provides an affordable, easy-to-use automation solution for low to medium throughput users of the Lonza PyroGene® rFC Endotoxin Detection Assay.
- High quality, reproducible results can be achieved with the ASSIST PLUS, eliminating the risk of costly and time-consuming retests.
- The ease and flexibility of VIALAB software allows users to customize plate layouts or set up partial plates using the same labware defined in this application note.

Materials

Manufacturer	Part Number	Description	Link
INTEGRA Biosciences	4505	ASSIST PLUS base unit	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus
INTEGRA Biosciences	4531	D-ONE Single channel 0.5-300 µl electronic pipetting module	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus#parts-and-numbers
INTEGRA Biosciences	4723	VOYAGER 8 channel 300 µl electronic pipette	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus#parts-and-numbers
INTEGRA Biosciences	4535	Tip deck for D-ONE on ASSIST PLUS	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus#parts-and-numbers
INTEGRA Biosciences	4552	6 x 8 Tube Rack	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus
INTEGRA Biosciences	6484, Specific Lots	300 µl LONG sterile (certified endotoxin free) GRIPTIPS®	https://www.integra-biosciences.com/global/en/pipette-tips/griptip-selector-guide
INTEGRA Biosciences	4371, Specific lots	10 ml Reagent Reservoir, Sterile, SureFlo anti-sealing array, polystyrene (certified endotoxin free)	https://www.integra-biosciences.com/global/en/reagent-reservoirs/multichannel-reagent-reservoirs
Lonza	50-658U, 50-658NV	PyroGene Recombinant Factor C Endotoxin Detection Assay	https://bioscience.lonza.com/lonza_bs/US/en/recombinant-factor-c-assay?gclid=EAlalQobChMI9NiJ4NqB_gIVS21vBB1AEQcTEAAYASAAEgKQJvD_BwE
Lonza	25-345S	PyroWave XM Fluorescence Reader	https://bioscience.lonza.com/lonza_bs/US/en/Endotoxin-Detection/p/00000000000213090/Pyro-Wave%E2%84%A2-XM-Fluorescence-Reader
Lonza	25-611	WinKQCL Software	https://bioscience.lonza.com/lonza_bs/US/en/winkqcl-endotoxin-detection-and-analysis-software?gclid=CjwKCA-jwolqhBhAGEiwArXT7K3QLa3cp0ejDVyHeG75DKONHmU-9v9UiDI1LyMJ9etKiJGJT2dCr9xoCrhoQAvD_BwE
Lonza	N207	Pyrogen-free Test Tubes	https://www.lonzabioscience.com.au/product/pyrogen-free-dilution-tubes-13x100mm-without-cap/
Lonza	W50-640	LAL Reagent Water	https://bioscience.lonza.com/lonza_bs/CH/en/Endotoxin-Detection/p/00000000000187028/LAL-Reagent-Water
Corning	3603	96 well black polystyrene microplate	https://ecatalog.corning.com/life-sciences/b2b/US/en/Microplates/Assay-Microplates/96-Well-Microplates/Corning%C2%AE-96-well-Black-Clear-and-White-Clear-Bottom-Polystyrene-Microplates/p/3603
IKA		Vortex 2	https://www.ika.com/en/Products-LabEq/Shakers-pg179/Vortex-2-25000258/

INTEGRA Biosciences AG
7205 Zizers, Switzerland
T +41 81 286 95 30
F +41 81 286 95 33
info@integra-biosciences.com

INTEGRA Biosciences Corp.
Hudson, NH 03051, USA
T +1 603 578 5800
F +1 603 577 5529
info-us@integra-biosciences.com

INTEGRA Biosciences Deutschland GmbH
35444 Biebertal, Deutschland
T +49 6409 81 999 15
F +49 6409 81 999 68
info-de@integra-biosciences.com

INTEGRA Biosciences SAS
95062 Cergy-Pontoise Cedex 1, France
T +33 (0)1 34 30 76 76
F +33 (0)1 34 30 76 79
info-fr@integra-biosciences.com

INTEGRA Biosciences Ltd.
Thatcham, Berks RG19 4EP, UK
T: +44 1635 797000
F: +44 1635 797001
info-uk@integra-biosciences.com

INTEGRA Biosciences Nordic ApS
Vallensbækvej 22A 3TV
Brøndby 2605, Denmark
T + 45 3173 5373
info-nordic@integra-biosciences.com

INTEGRA Biosciences (Shanghai) Co., Ltd.
中国上海自由贸易试验区环科路515号1110室
邮编: 201315
电话: +86 21 5844 7203
info-cn@integra-biosciences.com

インテグラ・バイオサイエンス(株)
〒101-0031
東京都千代田区東神田 1-5-6
東神田MK第五ビル 3階
T 03-5962-4936
F 03-5962-4937
info-jp@integra-biosciences.com